



## Protein Extraction from Cells Cultured in Alvetex® Scaffold 96-Well Plate Format

### Introduction

This protocol describes the use of a commercial protein extraction reagent to obtain protein samples from cells cultured in Alvetex Scaffold 96-well plate format (AVP009). Using this method, quantitative protein extraction is achievable over wide ranging cell densities. Resulting protein extracts are compatible with common downstream techniques, including Bradford and BCA protein assays, SDS-PAGE and western blotting, non-denatured immunological applications (e.g. immunoprecipitation), and enzymatic assays.

### Method

This is an example protocol, based on 3D culture of MDA-MB-231 breast cancer cells in Alvetex Scaffold in 96-well plates. Total protein was extracted using the Mammalian Protein Extraction Reagent (M-PER) from Pierce (product code 78503, [http://www.piercenet.com/browse.cfm?fldID=0601\\_0420](http://www.piercenet.com/browse.cfm?fldID=0601_0420)).

1. Withdraw culture medium from each well of the Alvetex Scaffold 96-well plate.
2. Rinse cells by carefully adding 200 DI PBS per well, aspirate to waste.
3. Add 100 µl M-PER reagent to each well and place plate on a shaking platform (100 rpm for 10 minutes).



**Note:** The M-PER reagent can be supplemented with protease inhibitors (e.g. Pierce, 87785) to minimise protein degradation upon release of cellular proteases.

4. Agitate the liquid in each well by pipetting up and down, and transfer to 0.5 ml microcentrifuge tubes.
5. Centrifuge samples at 14,000 x g for 5 minutes to remove cell debris.
6. Transfer supernatants to clean tubes and store at -80 °C.

### Example data

MDA-MB-231 breast carcinoma cells (ATCC, HTB-26) were routinely maintained in T-75 flasks. MDA-MB-231 culture media consisted of: Leibovitz's L-15 medium (ATCC, 30-2008) supplemented with 10 % v/v FBS and 100 U/ml Penicillin/Streptomycin. Cells were seeded into Alvetex Scaffold 96-well plates (AVP009) at seeding densities ranging from  $25 \times 10^3$  to  $200 \times 10^3$  cells in a total media volume of 200 µl per well. Plates were incubated for 24 hours at 37 °C with 0 % CO<sub>2</sub>, after which cultures were processed according to the protocol above.

# Protein Extraction from Cells Cultured in Alvetex® Scaffold 96-Well Plate Format

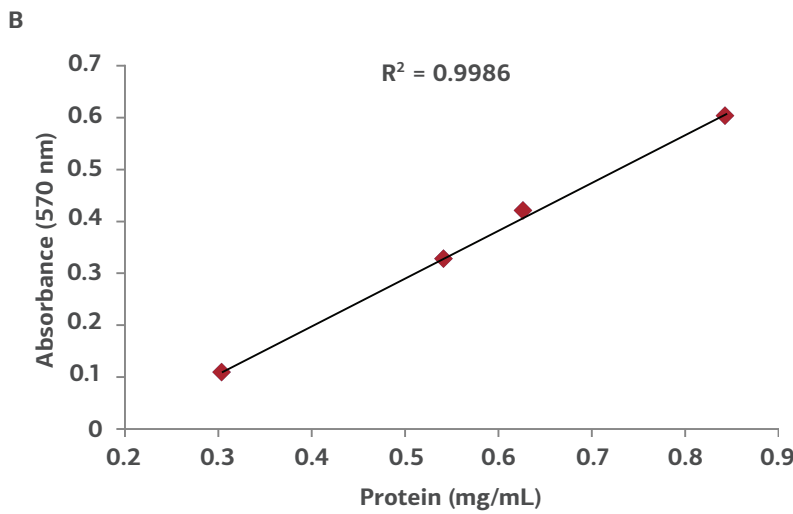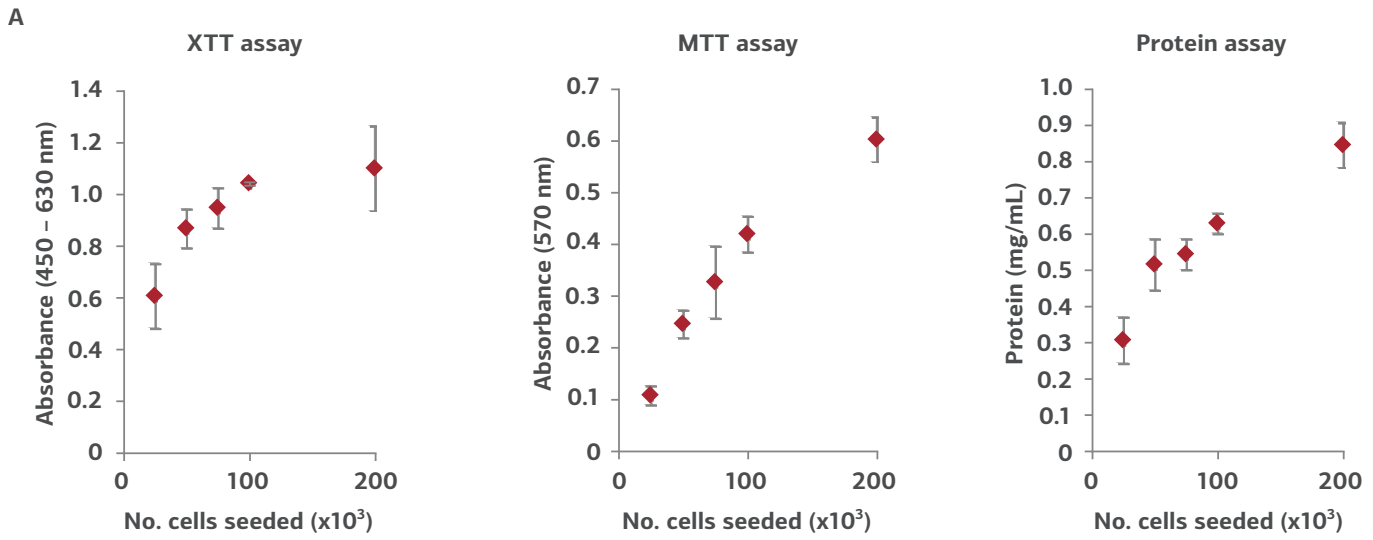

**C**

No. cells seeded per well	Protein yield per 3D culture (µg/ml)
25,000	30.5
50,000	51.4
75,000	54.3
100,000	62.8
200,000	84.4

**FIGURE 1.** Protein extraction from Alvetex Scaffold 96-well plate. MDA-MB-231 cells were seeded into Alvetex Scaffold 96-well plates at densities of 25, 50, 75, 100, and 200 x 10<sup>3</sup> cells in 200 µl per well. Cell viability was measured using two methods (XTT and MTT assays) for comparison with protein concentrations in extracts obtained via the protocol above (A) (n=3, mean ± SD). Pairwise plot of MTT viability and protein concentration (B) reveals a linear relationship between cell number and protein yield (R<sup>2</sup> = 0.9986). Protein yields ranged from approximately 30 µg/well for 25 x 10<sup>3</sup> cells seeded, to approximately 85 µg/well for 200 x 10<sup>3</sup> cells (C). Protein was estimated using the Protein Assay Reagent (Bio-Rad, 500-0006) and a BSA standard curve. For protocols for XTT and MTT viability assays in Alvetex Scaffold 96-well plates see: <http://reinnervate.com/alvetex/alvetex-documentation/protocols/>. Approximate protein yields in extracts from Alvetex Scaffold cultures of HepG2 hepatocytes, using BSA as a standard reference. Values shown represent the mean of triplicate readings per sample (±SD).

For research use only. Not for use in diagnostic procedures.



## Protein Extraction from Cells Cultured in Alvetex® Scaffold 96-Well Plate Format

### Technical Support

To obtain the latest technical support information, please visit [www.reinnervate.com](http://www.reinnervate.com). You can have 24-hour access to the following information:

- Access to telephone and fax information for order support
- Search through the Reinnervate protocol video gallery
- Search product spec sheets, protocols, applications notes and product support docs
- Obtain information on customer training and workshops
- Submit a question directly to Technical Support at [tech.support@reinnervate.com](mailto:tech.support@reinnervate.com)

---

### Ordering Information

For the latest products and services available through Reinnervate, please visit [www.reinnervate.com](http://www.reinnervate.com).

#### Reinnervate

NETPark Incubator, Thomas Wright Way,  
Sedgefield, Co. Durham, TS21 3FD  
UNITED KINGDOM

T: +44 (0) 1740 625 266

F: +44 (0) 1740 625 251

E: [info@reinnervate.com](mailto:info@reinnervate.com)

---

### Trademarks

Unless otherwise noted, all trademarks are the property of Reinnervate Ltd.  
©2015 Reinnervate Ltd. All rights reserved.