



AlamarBlue® Cell Viability Assay of upcyte® hepatocyte Cultures Grown on Alvetex® Scaffold in 3D

Introduction

The AlamarBlue® assay is based on the biological reduction of resazurin to the fluorescent molecule, resorufin. Resazurin is nontoxic and will readily penetrate into viable cells whereupon it is metabolised. The read-out options for the assay are absorbance or fluorescence, the latter being the more sensitive. The resazurin substrate is blue in colour and virtually non-fluorescent, resulting in very low background values in the fluorescent read out.

The amount of resorufin product increases with time and deviation from linearity can be anticipated at later time points. Therefore, it is recommended that for each cell type the linear relationship between both cell number and time, and signal produced should be first established in order to investigate the limits of the assay.

Methods

These are example protocols which may require further optimisation depending on the rate of cell growth of a particular cell type on Alvetex Scaffold. The protocols describe cell viability analysis of upcyte® hepatocytes (Medicyte, www.medicyte.com) using the AlamarBlue® Assay (Invitrogen, DAL1010, 1025, 1100). Upcyte® cells were thawed, immediately seeded onto the Alvetex Scaffolds and medium added after a 3 hour attachment period. At various time-points, the growth and viability of the upcyte® hepatocytes were monitored as described below.

Method 1: Performing the AlamarBlue® assay on dismantled Alvetex Scaffold discs

1. Prepare sufficient AlamarBlue® solution for the experiment, including controls, by diluting 1/10 in appropriate phenol red-free medium.
2. Remove the Alvetex Scaffold disc from the culture vessel and dismantle inserts where applicable. For 22 mm Alvetex Scaffold discs, transfer to the wells of a clean 12-well plate. For 15 mm Alvetex Scaffold discs, transfer to the wells of a clean 24-well plate.
3. For 22 mm Alvetex Scaffold discs add 0.8 ml of diluted AlamarBlue® solution to each well. Also add 0.8 ml diluted AlamarBlue® solution to empty control wells containing no cells. For 15 mm Alvetex Scaffold discs add 0.4 ml of diluted AlamarBlue® solution to each well. Also add 0.4 ml diluted AlamarBlue® solution to empty control wells containing no cells.
4. Incubate samples at 37 °C and 5 % CO₂ for 1 hour.
5. Transfer 100 µl aliquots of each sample to replicate wells of a black 96-well plate.
6. Measure fluorescence at 590 nm, using an excitation wavelength of 530-560 nm.

Method 2: Performing the AlamarBlue® assay on Alvetex Scaffold *in situ* in well inserts

1. Prepare sufficient AlamarBlue® solution for the experiment, including controls, by diluting 1/10 in appropriate phenol red-free medium.



AlamarBlue® Cell Viability Assay of upcyte® hepatocyte Cultures Grown on Alvetex® Scaffold in 3D

2. Remove the Alvetex Scaffold insert from the culture vessel and transfer to a clean plate. For 22 mm Alvetex Scaffold inserts, transfer to a clean 6-well plate. For 15 mm Alvetex Scaffold inserts, transfer to a clean 12-well plate.
3. For 22 mm Alvetex Scaffold inserts add 7 ml of diluted alamarBlue® solution to each well. For 15 mm Alvetex Scaffold inserts add 3 ml of diluted alamarBlue® solution to each well. Add the same corresponding volume of diluted alamarBlue® solution to empty control wells containing no cells.
4. Incubate samples at 37 °C and 5 % CO₂ for 3 hours.
5. Transfer 100 µl aliquots of each sample to replicate wells of a black 96-well plate.
6. Measure fluorescence at 590 nm, using an excitation wavelength of 530-560 nm.

Example Data Set

The example data shown below was obtained from cultures of upcyte® hepatocytes (Medicyte) grown in 15 mm Alvetex Scaffold presented in 12-well insert format.

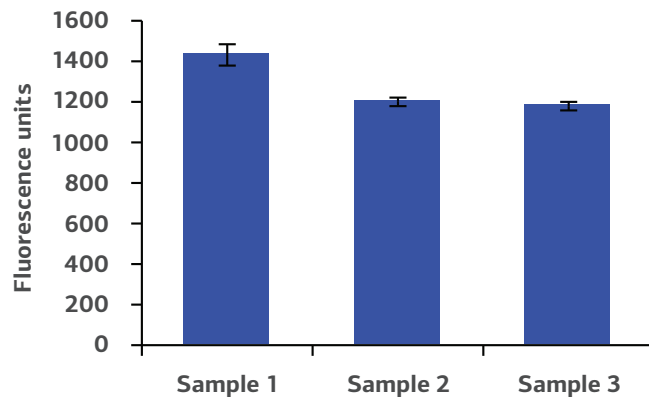


FIGURE 1. Biochemical analysis of cell viability using the alamarBlue® assay according to Method 2, above. Data from 3 sample replicates of upcyte® hepatocytes cultured in Alvetex Scaffold in 12-well insert format are shown (n=3, mean ± SD). Note good reproducibility between samples.



AlamarBlue® Cell Viability Assay of upcyte® hepatocyte Cultures Grown on Alvetex® Scaffold in 3D

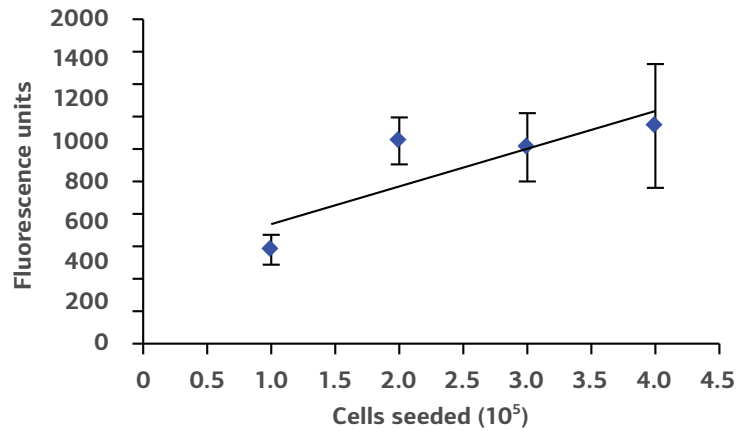


FIGURE 2. alamarBlue® assay of upcyte® hepatocytes cultured for 5 hours in Alvetex Scaffold discs presented in 12-well insert format (n=3, mean ± SD). Correlation is shown between cell number and fluorescent signal at seeding densities of 1.0, 2.0, 3.0 and 4.0 x 10⁵ cells per disc (R² > 0.71).

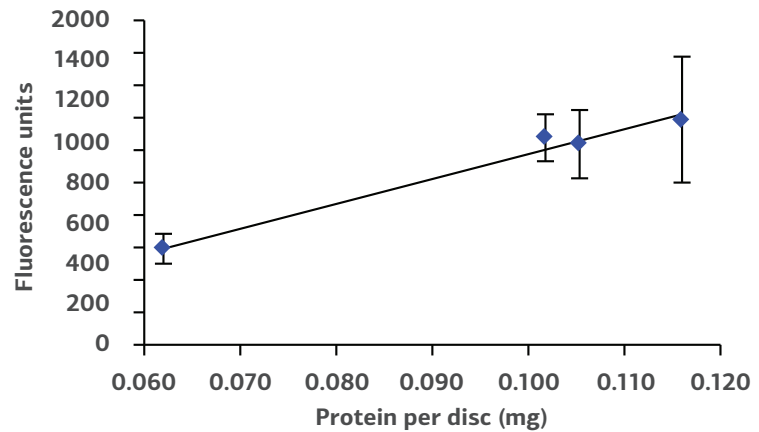


FIGURE 3. alamarBlue® assay of upcyte® hepatocytes cultured for 5 hours in Alvetex Scaffold discs presented in 12-well insert format (n=3, mean ± SD). Data normalised to protein extracted from Alvetex Scaffold discs seeded at 1.0, 2.0, 3.0 and 4.0 x 10⁵ cells. Good linearity was observed (R² = 0.98) between cellular protein and alamarBlue® fluorescence reading.

For research use only. Not for use in diagnostic procedures.



AlamarBlue® Cell Viability Assay of upcyte® hepatocyte Cultures Grown on Alvetex® Scaffold in 3D

Technical Support

To obtain the latest technical support information, please visit www.reinnervate.com. You can have 24-hour access to the following information:

- Access to telephone and fax information for order support
- Search through the Reinnervate protocol video gallery
- Search product spec sheets, protocols, applications notes and product support docs
- Obtain information on customer training and workshops
- Submit a question directly to Technical Support at tech.support@reinnervate.com

Ordering Information

For the latest products and services available through Reinnervate, please visit www.reinnervate.com.

Reinnervate

NETPark Incubator, Thomas Wright Way,
Sedgefield, Co. Durham, TS21 3FD
UNITED KINGDOM

T: +44 (0) 1740 625 266

F: +44 (0) 1740 625 251

E: info@reinnervate.com

Trademarks

Unless otherwise noted, all trademarks are the property of Reinnervate Ltd.
©2015 Reinnervate Ltd. All rights reserved.