



Example Protocol for the Culture of the MG63 Cell Line on Alvetex® Scaffold in Well Insert Formats

Introduction

Alvetex Scaffold is currently available in four different cell culture formats: 24-well plate (AVP006), 12-well plate (AVP002), 6-well insert (AVP004), and 12-well insert (AVP005). 24-well and 12-well plates are suitable for shorter term cultures and for applications where limited cell penetration into the scaffold is required. Well insert formats generally support longer term cultures and deeper cell penetration into the scaffold. They also provide for conveniently tailored media set ups (see Quick Start Protocol). 6-well inserts can be placed in conventional 6-well plates, while 12-well inserts can be placed in either 6-well plates or 12-well plates, depending on media requirements. Alternatively, both insert types can be housed in the dedicated Well Insert Holder in Deep Well Petri Dish (AVP015) to allow for increased media volumes and prolonged cell culture. The availability of two different Alvetex Scaffold sizes enables choice on the basis of desired culture size and cell expenditure.

Methods

Preparation for 3D Cell Culture on Alvetex Scaffold

1. MG63 cells were routinely maintained in T-75 flasks.

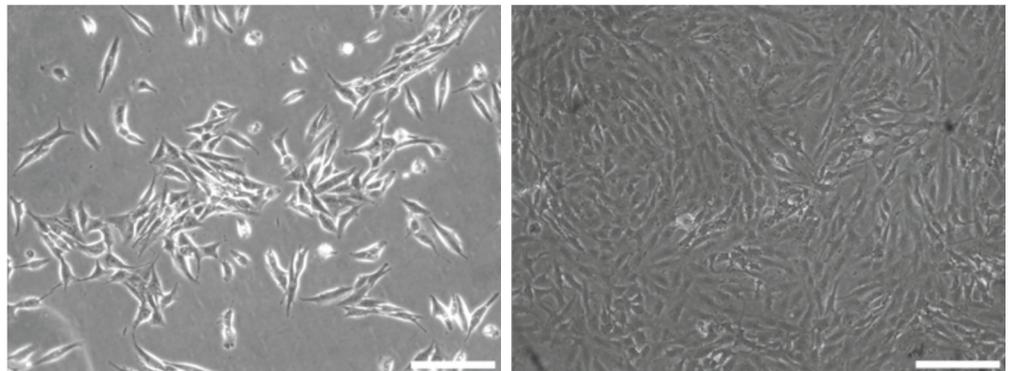


FIGURE 1. Phase contrast micrographs of MG63 cells grown in conventional 2D culture plates. Images show cells at low (left) and high (right) confluency. Scale bars: 100 µm.

2. Complete growth media consisted of: DMEM High glucose supplemented with 10 % v/v heat inactivated FBS, 1X non-essential amino acids, 2 µm L-glutamine and 100 U/ml Penicillin/ Streptomycin.
3. Cells were harvested by trypsinisation and centrifuged for 5 minutes (1000 rpm). The supernatant was discarded and the cell pellet was re-suspended in an appropriate volume of media for cell counting by Trypan Blue.
4. Cells were re-suspended at a concentration of 1.0×10^7 cells/ml for seeding.



Example Protocol for the Culture of the MG63 Cell Line on Alvetex® Scaffold in Well Insert Formats

12-well Insert Format, AVP005

1. Alvetex Scaffold 12-well inserts in 6-well plate format were prepared for seeding by dipping in 70 % ethanol and washed twice with media (7 ml per well).
2. 50 µl of the cell suspension was added to the centre of the Alvetex Scaffold disc, which was equivalent to 0.5×10^6 cells per well.
3. 5 ml of media was added to each well taking care not to allow media over the windows of the insert, i.e. independent feeding from below. The plate was incubated overnight at 37 °C with 5 % CO₂ to allow the cells to settle into the scaffold.
4. A further 5 ml of media was added to each well the following morning taking care not to dislodge cells from Alvetex Scaffold.
5. Plates were re-incubated and maintained by complete media exchange after every 2-3 days.

Note: This method can be applied to the use of Alvetex Scaffold in 6-well insert format, AVP004. Adjust cell seeding and media volumes according to the guidelines provided in the Quick Start Protocol.

Example Data: 12-well insert format, AVP005

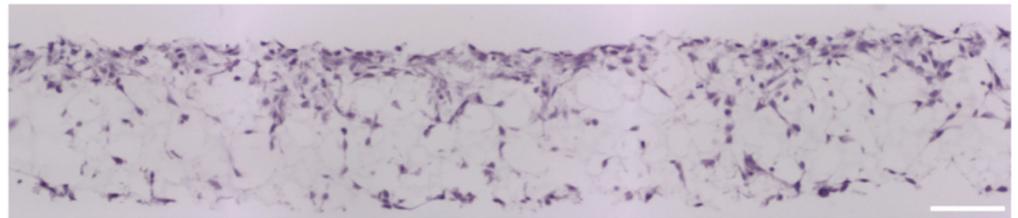


FIGURE 2. Brightfield micrographs showing the structure of MG63 cells cultured for 7 days on 15 mm diameter Alvetex Scaffold discs presented in 12-well insert in 6-well plate format. Cells were fixed, embedded in paraffin wax, sectioned (10 µm) and counterstained with haematoxylin and eosin. Scale bar 100 µm.



Example Protocol for the Culture of the MG63 Cell Line on Alvetex® Scaffold in Well Insert Formats

Technical Support

To obtain the latest technical support information, please visit www.reinnervate.com. You can have 24-hour access to the following information:

- Access to telephone and fax information for order support
- Search through the Reinnervate protocol video gallery
- Search product spec sheets, protocols, applications notes and product support docs
- Obtain information on customer training and workshops
- Submit a question directly to Technical Support at tech.support@reinnervate.com

Ordering Information

For the latest products and services available through Reinnervate, please visit www.reinnervate.com.

Reinnervate

NETPark Incubator, Thomas Wright Way,
Sedgefield, Co. Durham, TS21 3FD
UNITED KINGDOM

T: +44 (0) 1740 625 266

F: +44 (0) 1740 625 251

E: info@reinnervate.com

Trademarks

Unless otherwise noted, all trademarks are the property of Reinnervate Ltd.
©2015 Reinnervate Ltd. All rights reserved.