



Example protocol for the Culture of the 3T3 Cell Line on Alvetex® Scaffold in Well Plate and Well Insert Formats

Introduction

Alvetex Scaffold is currently available in four different cell culture formats: 24-well plate (AVP006), 12-well plate (AVP002), 6-well insert (AVP004), and 12-well insert (AVP005). 24-well and 12-well plates are suitable for shorter term cultures and for applications where limited cell penetration into the scaffold is required. Well insert formats generally support longer term cultures and deeper cell penetration into the scaffold. They also provide for conveniently tailored media set ups (see Quick Start Protocol). The availability of two different well insert formats enables choice on the basis of desired culture size and cell expenditure. 6-well inserts can be placed in conventional 6-well plates, while 12-well inserts can be placed in either 6-well plates or 12-well plates, depending on media requirements. Alternatively, both insert types can be housed in the dedicated Well Insert Holder in Deep Well Petri Dish (AVP015) to allow for increased media volumes and prolonged cell culture.

Methods

Preparation for 3D cell culture on Alvetex Scaffold

1. 3T3 cells were routinely maintained in T-75 flasks.

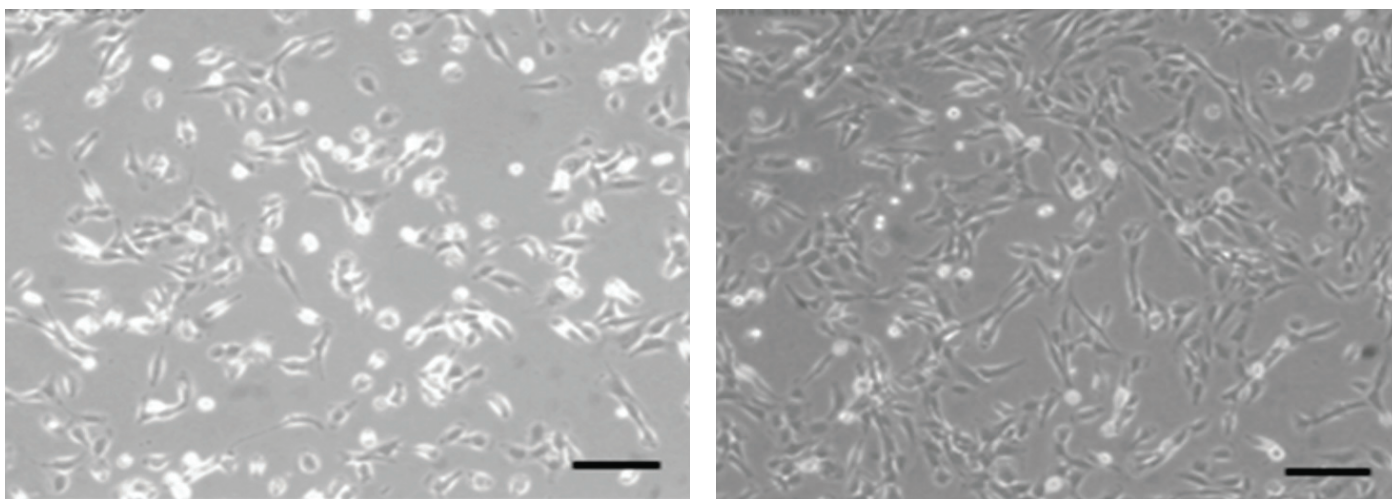


FIGURE 1. Phase contrast micrographs of 3T3 cells grown in conventional 2D culture plates. Images show cells at low (left) and high (right) confluency. Scale bars: 200 µm.

2. Complete media consisted of: Dulbecco's Modified Eagles Medium (DMEM) supplemented with 10% v/v FBS, 2 mM L-glutamine and 100 U/ml Penicillin/Streptomycin.
3. Cells were harvested by trypsinisation and centrifuged for 5 minutes (1000 rpm). The supernatant was discarded and the cell pellet was re-suspended in an appropriate volume of media for cell counting by Trypan Blue.
4. Cells were re-suspended at a concentration of 8×10^6 cells/ml for seeding.



Example protocol for the Culture of the 3T3 Cell Line on Alvetex® Scaffold in Well Plate and Well Insert Formats

5. Follow one of the below methods according to your choice of Alvetex Scaffold format.

12-well Plate Format, AVP002

1. Alvetex Scaffold 12-well plates were prepared for seeding with 70% ethanol (2 ml per well) and media washes (twice with 3 ml of media each).
2. 125 µl of the cell suspension was added to the centre of the Alvetex Scaffold disc, which was equivalent to 1×10^6 cells per well.
3. The plate was incubated 3 hours at 37 °C with 5% CO₂ to allow the cells to settle into the scaffold.
4. 4 ml of media was added to each well taking care not to dislodge cells from Alvetex Scaffold.
5. Plates were re-incubated and maintained by complete media exchange after every 1-2 days.

Note: This method can be applied to the use of Alvetex Scaffold in 24-well plate format, AVP006. Adjust cell seeding and media volumes according to the guidelines provided in the Quick Start Protocol.

6-well Insert Format, AVP004

1. Alvetex Scaffold 6-well inserts were prepared for seeding by dipping in 70 % ethanol and washed twice with media (7 ml per well).
2. 125 µl of the cell suspension was added to the centre of the Alvetex Scaffold disc, which was equivalent to 1×10^6 cells per well.
3. The plate was incubated for 60 minutes at 37 °C with 5 % CO₂ to allow the cells to settle into the scaffold.
4. 10 ml of media was added to each well taking care not to dislodge cells from Alvetex Scaffold.
5. Plates were re-incubated and maintained by complete media exchange after every 2-3 days.

Note: This method can be applied to the use of Alvetex Scaffold in 12-well insert format, AVP005. Adjust cell seeding and media volumes according to the guidelines provided in the Quick Start Protocol.

Example protocol for the Culture of the 3T3 Cell Line on Alvetex® Scaffold in Well Plate and Well Insert Formats



Example Data: 12-well plate format, AVP002

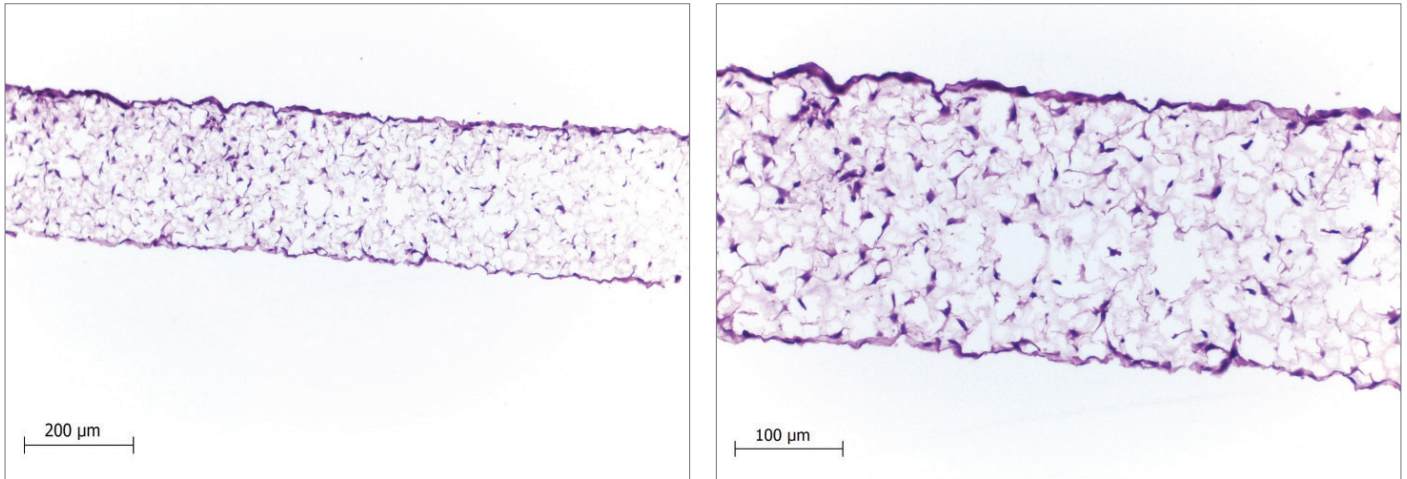


FIGURE 2. Brightfield micrographs showing the structure of 3T3 cells cultured for 7 days on 22 mm diameter Alvetex Scaffold discs presented in 2-well plate format. Cells were fixed, embedded in paraffin wax, sectioned (10 µm) and counterstained with haematoxylin and eosin.

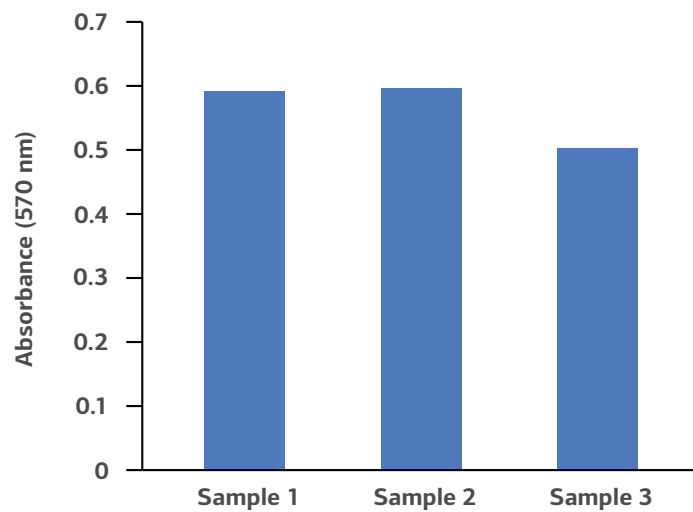


FIGURE 3. Biochemical analysis of cell viability using a standard MTT assay. Data from 3 sample replicates of 3T3 cells are shown. Each well was sampled in duplicate with mean value shown. Cells were cultured for 3 days on 22 mm Alvetex Scaffold discs presented in the 12-well plate format.

For research use only. Not for use in diagnostic procedures.

Example protocol for the Culture of the 3T3 Cell Line on Alvetex® Scaffold in Well Plate and Well Insert Formats



Example Data: 12-well plate format, AVP002

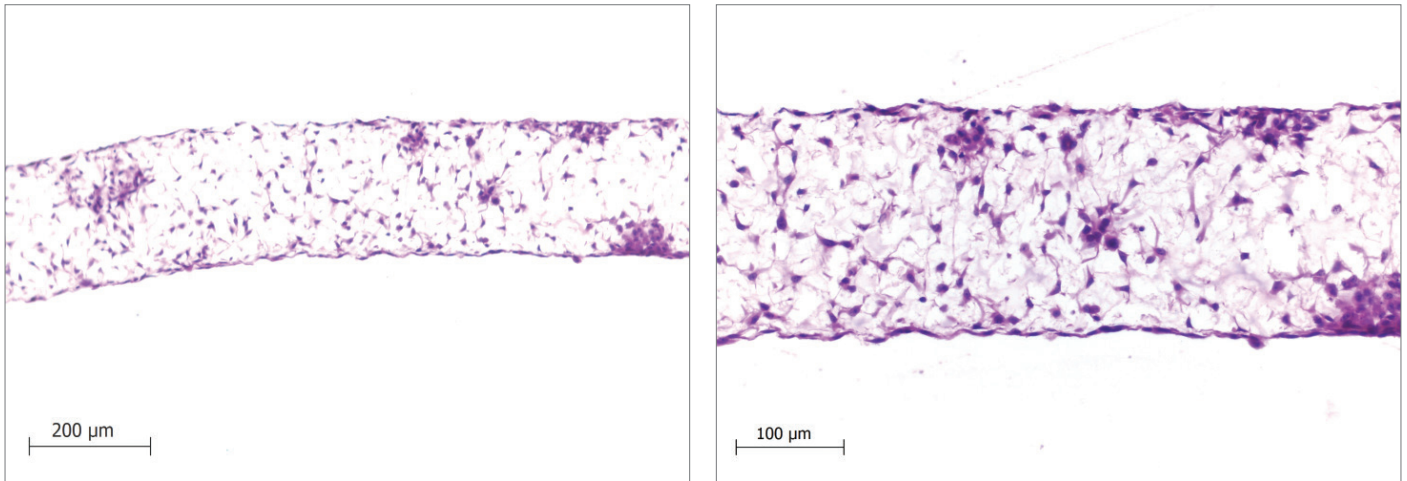


FIGURE 4. Brightfield micrographs showing the structure of 3T3 cells cultured for 7 days on 22 mm diameter Alvetex Scaffold discs presented in 6-well insert in 6-well plate format. Cells were fixed, embedded in paraffin wax, sectioned (10 µm) and counterstained with haematoxylin and eosin.

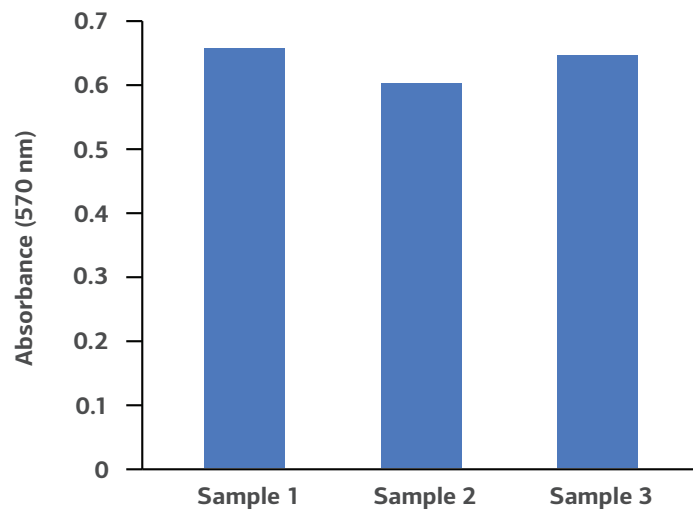


FIGURE 5. Biochemical analysis of cell viability using a standard MTT assay. Data from 3 sample replicates of 3T3 cells are shown. Each well was sampled in duplicate with mean value shown. Cells were cultured for 3 days on 22 mm Alvetex Scaffold discs presented in the 6-well inserts in 6-well plate format.

For research use only. Not for use in diagnostic procedures.



Example protocol for the Culture of the 3T3 Cell Line on Alvetex® Scaffold in Well Plate and Well Insert Formats

Technical Support

To obtain the latest technical support information, please visit www.reinnervate.com. You can have 24-hour access to the following information:

- Access to telephone and fax information for order support
- Search through the Reinnervate protocol video gallery
- Search product spec sheets, protocols, applications notes and product support docs
- Obtain information on customer training and workshops
- Submit a question directly to Technical Support at tech.support@reinnervate.com

Ordering Information

For the latest products and services available through Reinnervate, please visit www.reinnervate.com.

Reinnervate

NETPark Incubator, Thomas Wright Way,
Sedgefield, Co. Durham, TS21 3FD
UNITED KINGDOM

T: +44 (0) 1740 625 266

F: +44 (0) 1740 625 251

E: info@reinnervate.com

Trademarks

Unless otherwise noted, all trademarks are the property of Reinnervate Ltd.
©2015 Reinnervate Ltd. All rights reserved.