


Xeno-free, serum-free culture medium for human iPS and ES cells

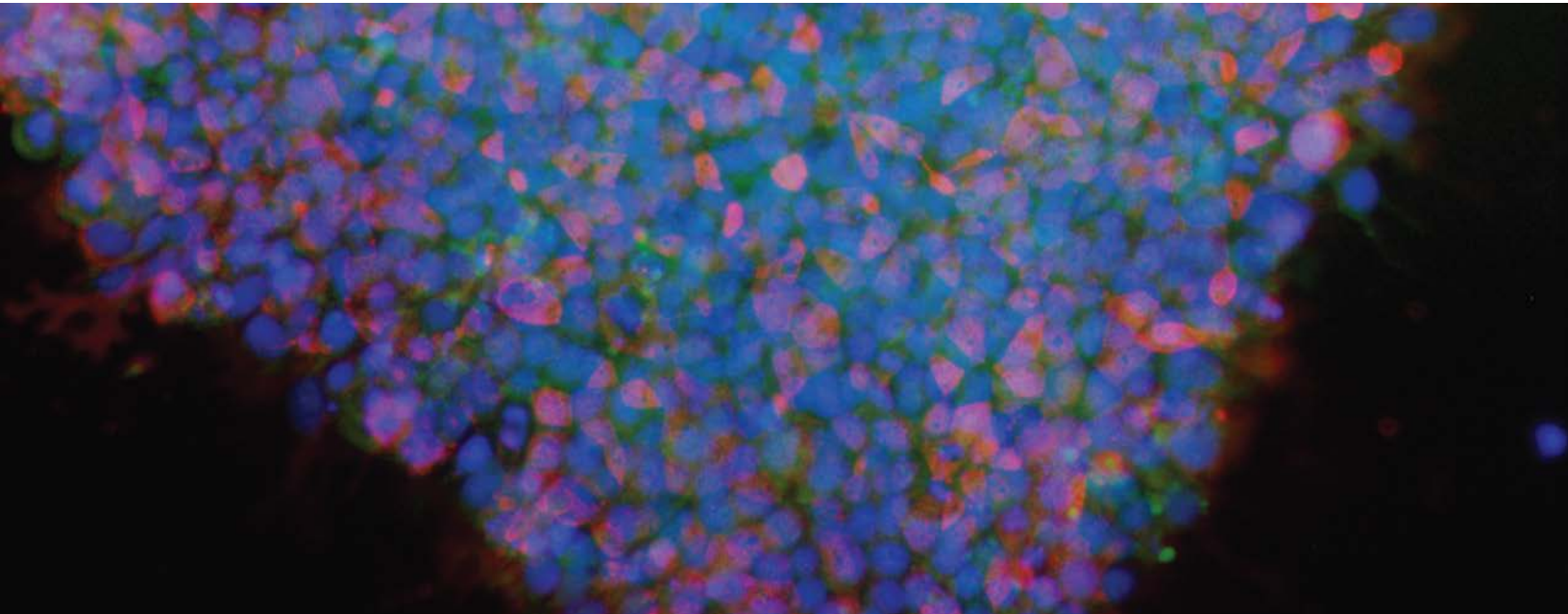

Expanding the culture of excellence

- **Defined, serum-free, xeno-free**
- **Complete, ready-to-use**
- **Superior proliferation of undifferentiated hES and hiPS cells**
- **Stable pluripotency and genotype over long-term culture**
- **Extensively tested and widely referenced**

NutriStem[®] XF/FF Culture Medium is a defined, xeno-free, serum-free medium designed to support the growth and expansion of human induced pluripotent stem (hiPS) and human embryonic stem (hES) cells in a feeder-free environment. NutriStem XF/FF Culture Medium offers the ability to culture human pluripotent cells without the need for high levels of bFGF and other stimulatory growth factors or cytokines. The low-protein formulation contains only the most essential components required for maintenance of hES and hiPS cells, providing a simplified medium and maintaining the cells' full differentiation potential.

The defined, xeno-free formulation of NutriStem XF/FF Culture Medium provides consistent media performance and predictable cellular behavior, as well as increased reproducibility in long-term culture (over 50 passages). In addition, cells cultured in NutriStem XF/FF Culture Medium show superior attachment and proliferation rates, making this medium ideal for high-throughput screening applications.

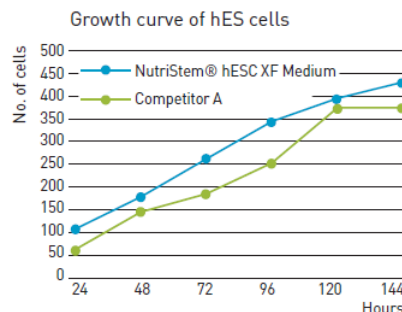


Figure 1: NutriStem[®] hPSC XF Medium enables excellent proliferation of undifferentiated hES and hiPS cells. Proliferation of H1 hES cells cultured in Corning Matrigel-coated 96-well plates in NutriStem XF/FF Medium and the leading competing medium for feeder-free culture. Medium was changed and proliferation was assessed every 24 hours in culture.

Normal cell morphology and functional assessment of pluripotency

The formation of compact colonies of cells with a high nucleus-to-cytoplasm ratio, prominent nucleoli, and distinct colony borders are characteristic morphology traits of healthy undifferentiated hES and hiPS cells, and can be observed through a phase-contrast microscope (Figure 2). Human pluripotent stem cells hold the potential to differentiate into cell types of all three germ layers, i.e., endoderm, mesoderm, and ectoderm. This differentiation potential is assessed by the spontaneous differentiation within embryoid bodies cultured *in vitro* (Figure 3) and teratomas formed *in vivo* (Figure 4).

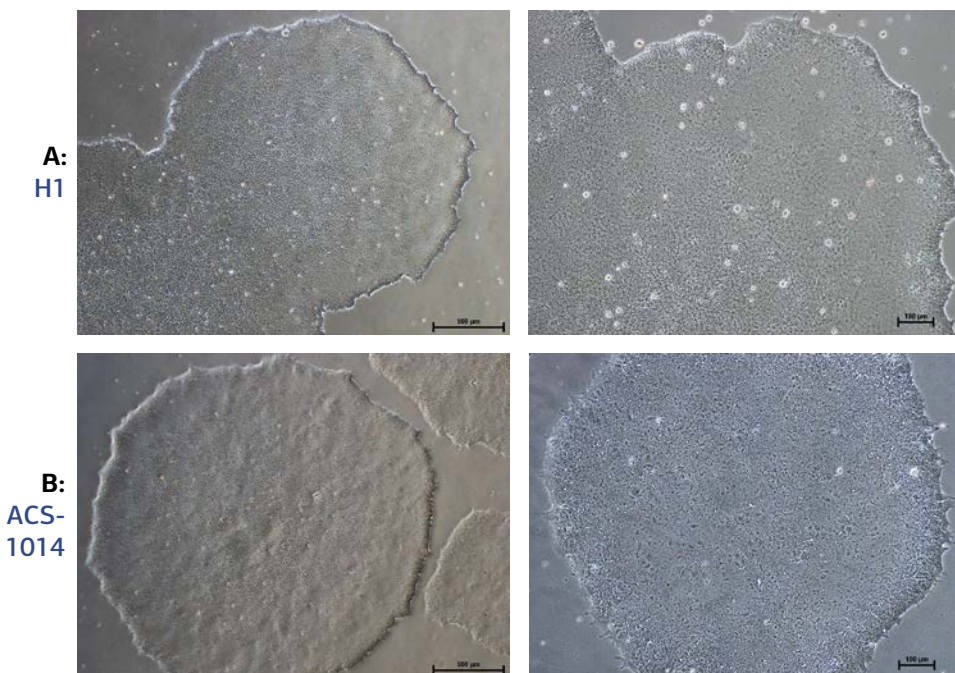


Figure 2: Normal Colony Morphology. H1 hES cells (panel A) and ACS-1014 hiPS cells (panel B) cultured in NutriStem XF/FF Medium on Corning Matrigel-coated plates display colony morphologies typical of normal feeder-free hES and hiPS cell cultures, including a uniform colony of tightly compacted cells and distinct colony edges.

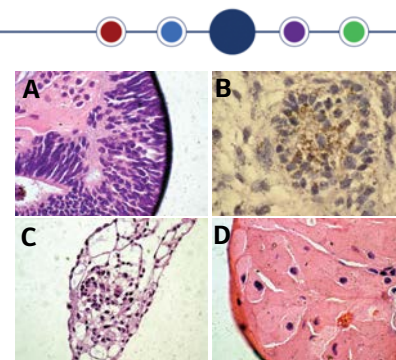


Figure 3: Embryoid Body Formation. Embryoid bodies (EBs) were generated from H9.2 hES cells cultured for 16 passages in NutriStem XF/FF Medium on Corning Matrigel matrix as an evaluation of pluripotency. The pluripotent H9.2 cells were suspended in serum-supplemented medium, where they spontaneously formed EBs containing cells of embryonic germ layers. The following cell types were identified by examination of the histological sections of 14-day-old EBs stained with H&E: (A) neural rosette (ectoderm), (B) neural rosette stained for tubulin, (C) primitive blood vessels (mesoderm), and (D) megakaryocytes (mesoderm).

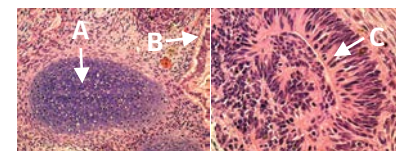


Figure 4: Teratoma Formation. H9.2 hES cells were cultured for 11 passages in NutriStem XF/FF Medium using a human foreskin fibroblast (HFF) feeder layer. The hES cells were subsequently injected into the hind leg muscle of SCID-beige mice for *in vitro* evaluation of pluripotency. The following tissues from all three germ layers were identified in H&E-stained histological sections of the teratoma 12 weeks post-injection: (A) cartilage (mesoderm), (B) epithelium (endoderm), and (C) neural rosette (ectoderm).

Ordering Information

Product Name	Cat. No.	Size
NutriStem XF/FF Culture Medium (BI Cat. No. 05-100-1)	01-0005	500 mL 100 mL
Related Product		
CryoStem™ Freezing Medium (BI Cat. No. 05-710-1)	01-0013	50 mL



AMERICAS • HUMAN TISSUE
9000 Virginia Manor Road, Suite 207
Beltsville, MD 20705, USA
Tel: (301) 470-3362
Email: info@bioserve.com
www.bioserve.com



AMERICAS • REAGENTS
4 Hartwell Place
Lexington, MA 02421, USA
Tel: (617) 245-0000
Email: info@stemgent.com
www.stemgent.com



ASIA PACIFIC
KDX Shin-yokohama381 Bldg9F
3-8-11, Shin-yokohama, Kohoku-ku,
Yokohama, Kanagawa 222-0033, Japan
Tel: +81 45 475 3887
Email: info_jp@reprocell.com
www.reprocell.com/en



EUROPE
NETPark Incubator
Thomas Wright Way
Sedgefield, Co. Durham
TS21 3FD, UNITED KINGDOM
Tel: +44 (0)1740 625 266
Email: info@reinnervate.com
www.reinnervate.com



**AMERICAS & EUROPE
HUMAN TISSUE ASSAYS**
Weipers Centre, Garscube Estate
Bearsden Road, Glasgow G61 1QH
United Kingdom
Tel: +44 (0)141 330 3831
Email: info@biopta.com
www.biopta.com