


Xeno-free, serum-free culture medium for human iPS and ES cells

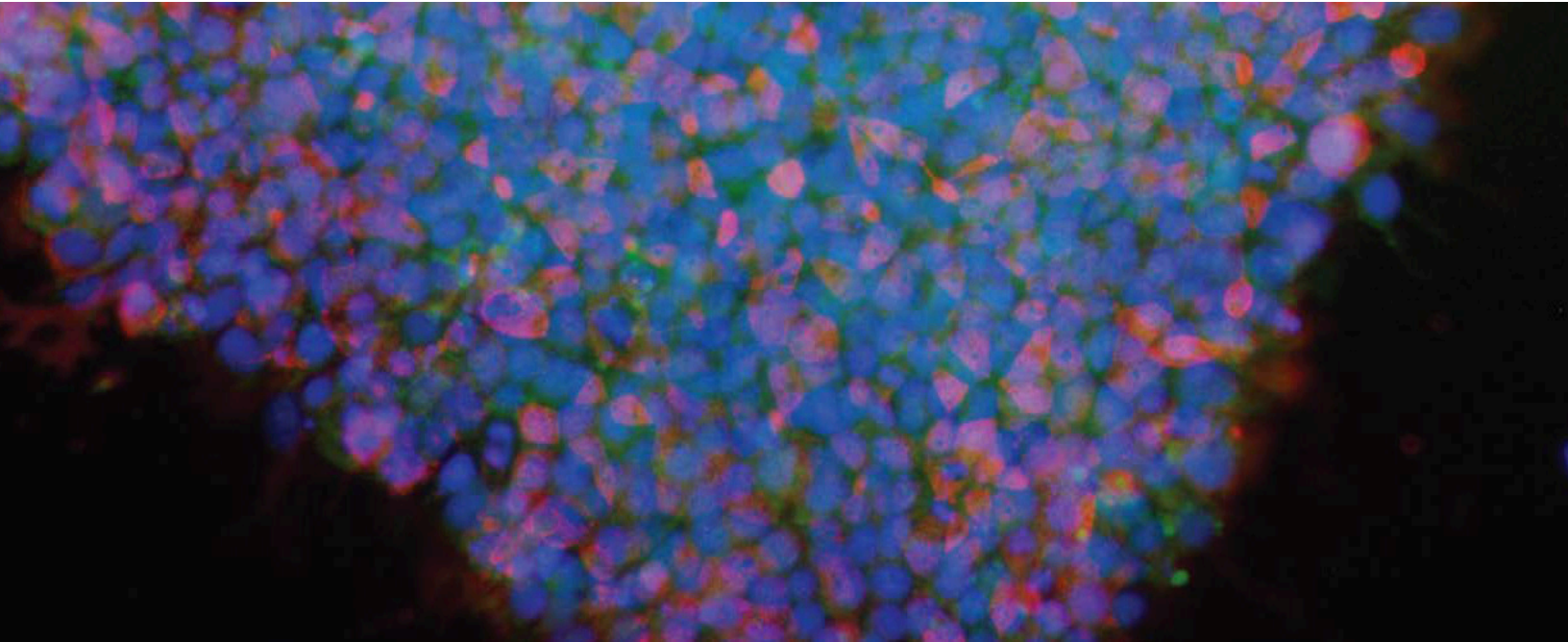

Expanding the culture of excellence

Biological Industries® NutriStem™ hPSC XF Culture Medium is a defined, xeno-free, serum-free medium designed to support the growth and expansion of human induced pluripotent stem (hiPS) and human embryonic stem (hES) cells in a feeder-free environment. NutriStem hPSC XF Culture Medium offers the ability to culture human pluripotent cells without the need for high levels of bFGF and other stimulatory growth factors or cytokines. The low-protein formulation contains only the most essential components required for maintenance of hES and hiPS cells, providing a simplified medium and maintaining the cells' full differentiation potential.

The defined, xeno-free formulation of NutriStem hPSC XF Culture Medium provides consistent media performance and predictable cellular behavior, as well as increased reproducibility in long-term culture (over 50 passages). In addition, cells cultured in NutriStem hPSC XF Culture Medium show superior attachment and proliferation rates, making this medium ideal for high-throughput screening applications.

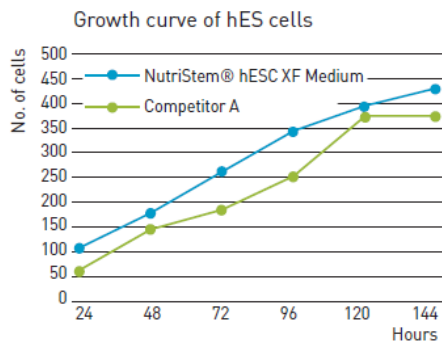


Figure 1. NutriStem® hPSC XF Medium enables excellent proliferation of undifferentiated hES and hiPS cells. Proliferation of H1 hES cells cultured in Corning Matrigel-coated 96-well plates in NutriStem hPSC XF Medium and the leading competing medium for feeder-free culture. Medium was changed and proliferation was assessed every 24 hours in culture.



- **Defined, serum-free, xeno-free**
- **Complete, ready-to-use**
- **Superior proliferation of undifferentiated hES and hiPS cells**
- **Stable pluripotency and genotype over long-term culture**
- **Extensively tested and widely referenced**

Biological Industries company name and logo and NutriStem and NutriFreez brand name are the property of Biological Industries Ltd., Israel. Company names, brands, logos and trademarks for REPROCELL, BioServe, Stemgent, Alvetex and Biopta are all the property of REPROCELL Inc. © 2019 REPROCELL, Inc. All rights reserved.

Normal cell morphology and functional assessment of pluripotency

The formation of compact colonies of cells with a high nucleus-to-cytoplasm ratio, prominent nucleoli, and distinct colony borders are characteristic morphology traits of healthy undifferentiated hES and hiPS cells, and can be observed through a phase-contrast microscope (**Figure 2**). Human pluripotent stem cells hold the potential to differentiate into cell types of all three germ layers, i.e., endoderm, mesoderm, and ectoderm. This differentiation potential is assessed by the spontaneous differentiation within embryoid bodies cultured *in vitro* (**Figure 3**) and teratomas formed *in vivo* (**Figure 4**).

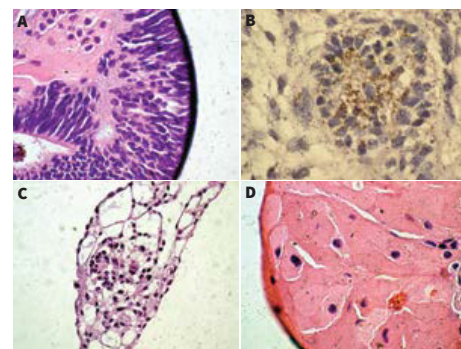


Figure 3. Embryoid Body Formation. Embryoid bodies (EBs) were generated from H9.2 hES cells cultured for 16 passages in NutriStem hPSC XF Medium on Corning Matrigel matrix as an evaluation of pluripotency. The pluripotent H9.2 cells were suspended in serum-supplemented medium, where they spontaneously formed EBs containing cells of embryonic germ layers. The following cell types were identified by examination of the histological sections of 14-day-old EBs stained with H&E: (A) neural rosette (ectoderm), (B) neural rosette stained for tubulin, (C) primitive blood vessels (mesoderm), and (D) megakaryocytes (mesoderm).

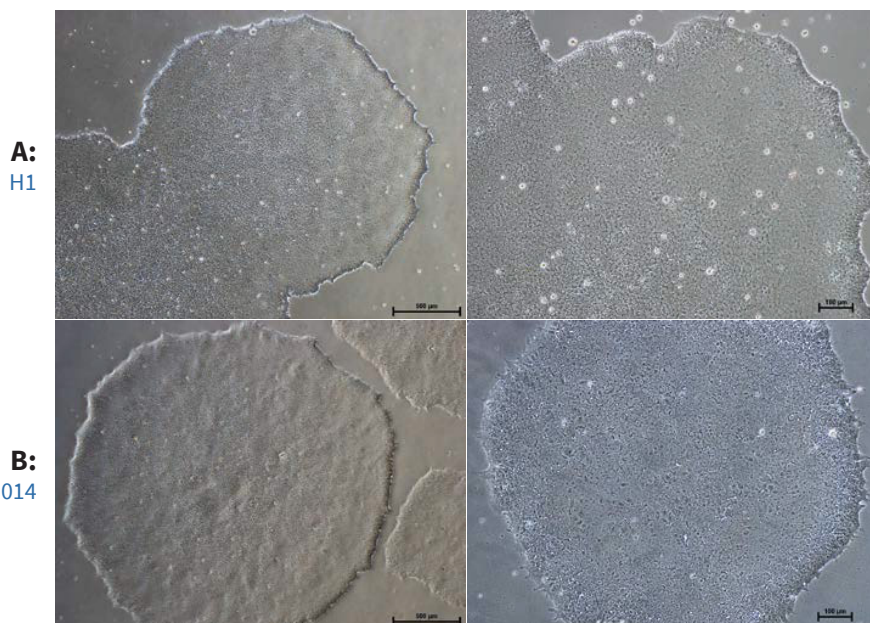


Figure 2. Normal Colony Morphology. H1 hES cells (panel A) and ACS-1014 hiPS cells (panel B) cultured in NutriStem hPSC XF Medium on Corning Matrigel-coated plates display colony morphologies typical of normal feeder-free hES and hiPS cell cultures, including a uniform colony of tightly compacted cells and distinct colony edges.

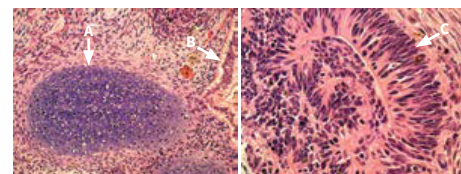


Figure 4. Teratoma Formation. H9.2 hES cells were cultured for 11 passages in NutriStem hPSC XF Medium using a human foreskin fibroblast (HFF) feeder layer. The hES cells were subsequently injected into the hind leg muscle of SCID-beige mice for *in vitro* evaluation of pluripotency. The following tissues from all three germ layers were identified in H&E-stained histological sections of the teratoma 12 weeks post-injection: (A) cartilage (mesoderm), (B) epithelium (endoderm), and (C) neural rosette (ectoderm).

Ordering Information

Product Name	Quantity	Cat. No.
NutriStem® XF hPSC Culture Medium (BI Cat. No. 05-100-1)	01-0005	500 mL
	01-0005-100	100 mL

Related Product

Product Name	Quantity	Cat. No.
NutriFreez™ D10 Cryopreservation Medium (BI Cat. No. 05-713-1E)	01-0020-50	50 mL

REPROCELL Europe Ltd

Thomson Pavillion, Todd Campus
West of Scotland Science Park
Acre Road
Glasgow, G20 0XA
UK

T: +44 (0)141 465 3460
E: info-europe@reprocell.com

REPROCELL USA Inc

9000 Virginia Manor Road
Suite 207
Beltsville, MD 20705
USA

T: +1 301 470 3362
E: info-us@reprocell.com

REPROCELL India Ltd

3-1-135/1A, CNR Complex
Mallapur
Hyderabad 500 076
Telangana
India

T: +91-40-27178178
E: Bhargavi.Gurram@reprocell.com

REPROCELL Inc (Japan)

MetLife Shin-yokohama 381, Bldg. 9F
3-8-11, Shin-yokohama
Kohoku-ku, Yokohama
Kanagawa 222-0033
Japan

T: +81 45 475 3887
E: info-asia@reprocell.com